

Photoresist Spinner Operation

- 1) Change the aluminum foil after each use. The build-up of photoresist can damage or impair proper operation.
- 2) Clean the spinner chuck with acetone after each use making sure to remove and subsequently replace the o-ring when cleaning.
- 3) Clean up any photoresist that has been spilled and wipe clean all surrounding surfaces.
- 4) Return the spinner chuck to its proper place when finished.
- 5) Report any problems or malfunction in this equipment immediately to the super-user and fill out an Incident Report located at the entrance of 1511.

- DISPOSE OF ANY PHOTORESIST CONTAMINATED MATERIALS IN THE PROPER WASTE CONTAINER. DO NOT THROW AWAY CONTAMINATED MATERIALS IN REGULAR TRASH.

Super user for Photoresist Spinner:

Aaron Birkbeck (alb@ucsd.edu)